
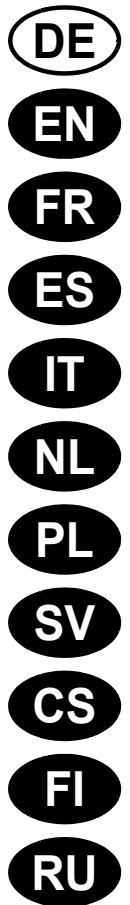
Installation and operating instructions
Notice de montage et de mise en service
Instrucciones de montaje y uso
Istruzioni per il montaggio e l'uso
Montage- en bedrijfsinstructies

Montaż i instrukcja obsługi
Monterings- og driftsvejledning
Návod pro montáž a provoz
Asennus- ja käyttöohje
Инструкция по монтажу и вводу в эксплуатацию

Best.-Nr.: 2000101103
FAID108

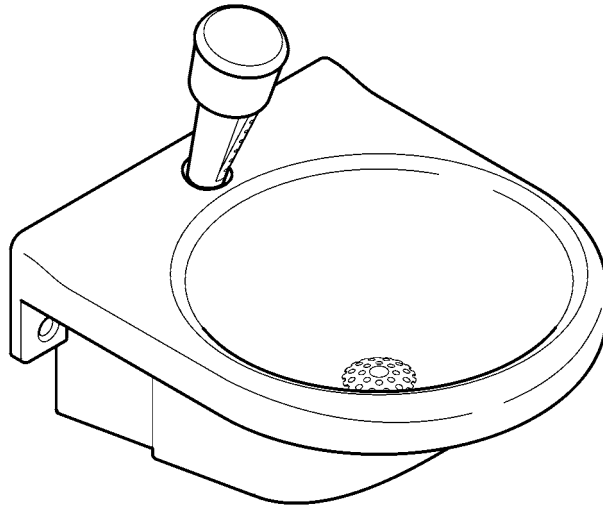EN

3

.....

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Eye and facial shower with water collection basin for wall-mounting

Please refer to the graphics in the German installation and operating instructions.

Table of contents

1. Abbreviations and units	3
2. Warranty	4
3. Important notes	4
Description of product	
4. Application	4
5. Technical specifications	4
6. Scope of delivery	4
7. Dimensions	4
8. Installation example	4
Assembly, function and commissioning	
9. Assembly with pressure reducer	5
10. Assembly without pressure reducer	5
11. Function	5
12. Adjust jet height	5
Maintenance	
13. Maintenance and care	6
14. Troubleshooting	6
15. Spare parts	6

1. Abbreviations and units

Best.-Nr.	Aquarotter order number
Conversion	1 mm = 0.03937 inches 1 inch = 25.4 mm

All length specifications in the graphics are in mm.

2. Warranty

Liability is accepted in accordance with the General Terms and Conditions of Business and Supply.

Use original spare parts only!

3. Important notes

- Assembly, commissioning and maintenance may only be carried out by a qualified technician in accordance with the instructions.
- The statutory regulations and conditions for technical connection to the local water supply companies must be adhered to.
- Prior to the assembly process, flush the pipework in accordance with DIN 1988.
- Mount the eye shower in accordance with EN 15154 Part 2.
- Replace the filter once a year.
- We reserve the right to make changes.

4. Application

Emergency showers are prescribed first aid installations for work stations where hazardous materials are being handled.

Eye and facial shower with water collection basin for wall mounting, including bracket for hand shower.

5. Technical specifications

Water connection:	G 3/8 B
Minimum flow pressure:	1.0 bar
Flow pressure range:	1-4 bar This range meets the requirements of EN 15154.
Volumetric flow:	Constant 6-9 l/s at a flow pressure of 1 bar

6. Scope of delivery

7. Dimensions

8. Installation example

9. Assembly with pressure reducer

For a flow pressure > 1.2 bar

9.1 Build the water drain.

9.2 Check the height of the water jet.

- ☞ If the jet height is outside the range of 100 - 300 mm, adjust the jet height (see [chapter 12.](#)).

10. Assembly without pressure reducer

For a flow pressure 1.0 - 1.2 bar

10.1 To mount the water collection basin, see [Work step 9.1 to 9.6.](#)

10.2 Mount the hose.

10.3 Build the water drain.

10.4 Check the height of the water jet.

- ☞ If the jet height is outside the range of 100 - 300 mm, adjust the jet height according to EN 15154 (see [chapter 12.](#)).

11. Function

11.1 Press the trigger button.

- The water will flow.

11.2 Press the red stop button.

- The water flow will stop.

- ☞ The fitting is not self-closing, so that the user can hold their eyes open with both hands!

- ☞ Depending on the length of the connection hose, the eye shower can also be used outside the water collection basin.

12. Adjust jet height

13. Maintenance and care

Check once a month that the eye shower is in working order.
(Guidelines for Laboratories BGI/GUV-I 850-0)

As a preventative measure to reduce microbiological contamination, it is recommended to replace the water in the fitting at shorter intervals by opening the valve.

Use suitable cleaning agents which are not harmful to the fitting in a proper manner and rinse with water after use. High-pressure cleaners must not be used for cleaning.

14. Troubleshooting

Fault	Cause	Remedy
Water does not flow	– Water supply interrupted – Dirty filter	⇒ Restore it ⇒ Clean and replace if necessary
Water volume too low	– Pressure reducer defective – Dirty filter	⇒ Replace it ⇒ Clean it
Jet height too high	– Jet height set incorrectly – Flow pressure too high	⇒ Set it ⇒ Assembly with pressure reducer
Jet height too low	– Supply pressure too low – Pressure reducer set incorrectly – Flow pressure too low	⇒ Check it ⇒ Set it ⇒ Assembly without pressure reducer

If you cannot rectify a fault or if it is not in the list, please contact our customer service department!

15. Spare parts

Bezeichnung	Best.-Nr.
1 Hand shower complete	2000105863
2 Pressure reducer	2000109525
3 Filter (packaging unit 10x)	2000109401
4 Dust cap with ball chain	2000105849
5 Spray head (packaging unit 4x)	2000104725
6 Non-return valve (packaging unit 10x)	2000104744
7 Seal (packaging unit 10x)	2000104668
8 Connection hose	2000105845
Maintenance set consisting of:	
O-rings	2000105864

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KWC ME LLC Ras Al Khaimah,
KWC DVS Ltd - Northern Office
KWC DVS Ltd - Southern Office
KWC Aquarotter GmbH
KWC Nordics Oy
KWC Austria GmbH